Aquatron

De-ionisers & Filters

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Assembly and Operating Instructions

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Aquatron De-ioniser ADH & Ion-exchange Cartridges ADI

Assembly Instructions

Your Aquatron De-ioniser is a highly versatile unit which can be assembled and operated in 3 modes:-

- 1 BENCH MOUNTED using the base support ring provided in the standard kit.
- 2 WALL MOUNTED using the optional wall bracket catalogue number AFD.
- 3 IN SERIES with other Aquatron De-ionisers and Filters using the optional plastic coupling catalogue number ALC.

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Bench Mounted Assembly

a Open the package containing the De-ioniser Housing ADH and identify the following components:

- De-ioniser lid
- Clear plastic bowl
- Base support ring
- Large spanner for bowl removal
- Small spanner for plumbing
- Plumbing components pack ADH/1
- Plumbing components pack ADH/2
- Plumbing components pack ADH/3

b Unscrew the clear plastic bowl from the de-ioniser lid using the large spanner.

NB: It is advisable to lubricate the rubber O ring on the lid with silicone grease. This permits easier assembly and disassembly.

- c Open the package containing the ion-exchange cartridges ADI. Select a cartridge and remove the plastic end caps. Place into bowl.
- d Refit the bowl and cartridge to the de-ioniser lid. Use the spanner to ensure a secure fit with no gap between the bowl and lid.
- e Firmly position the assembled unit into the base support ring.
- f Identify the "IN" and "OUT" ports on the De-ioniser lid.
- g Onto the "OUT" port assemble the plumbing components ADH/1 containing the water flow valve.
- h Onto the "IN" port assemble EITHER:
- the plumbing components ADH/2 containing a stainless flexible pipe and tap adapter.

OR - the plumbing components ADH/3 with plain olive connection.

NB: The choice may depend on the type and location of the mains water tap.

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
- i Use the small spanner provided to securely tighten all the components.
- j Now refer to OPERATING INSTRUCTIONS FOR AQUATRON DE-IONISER.

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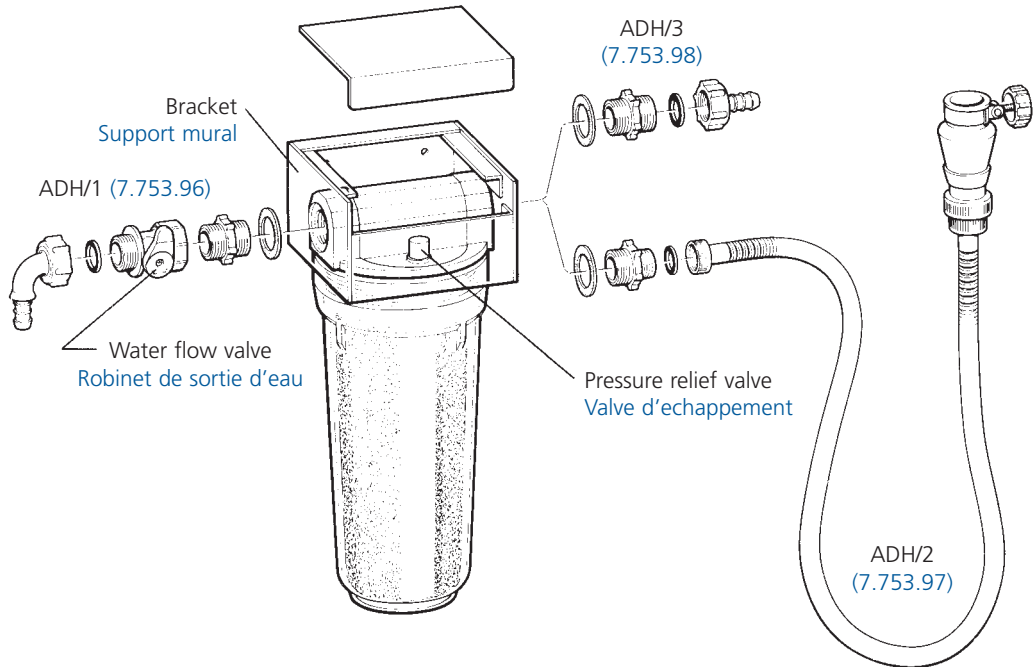Wall Mounted Assembly

For this mode of assembly you require:-

- Aquatron De-ioniser Housing ADH
 - Ion Exchange Cartridges ADI
 - Wall Mounting Bracket AFD
- a Fix the wall mounting bracket to a sound wall making sure the bracket is horizontal. Make sure the location of the bracket allows easy access for changing the ion-exchange cartridges.
 - b Unscrew the clear plastic bowl from the de-ioniser lid using the large spanner.
NB: It is advisable to lubricate the rubber O ring on the lid with silicone grease. This permits easier assembly and disassembly.
 - c Identify the "IN" and "OUT" ports on the de-ioniser lid.
 - d Position the lid inside the wall bracket with the ports in line with the bracket holes. Ensure the "OUT" port is facing the LEFT hole.
 - e Onto the "OUT" port assemble the plumbing components ADH/1 containing the water flow valve. Refer to Figure 2.
 - f Onto the "IN" port assemble EITHER:
 - the plumbing components ADH/2 containing a stainless flexible pipe and tap adapter.
 - OR - the plumbing components ADH/3 with plain olive connection.
NB the choice may depend on the type and location of the mains water tap.
 - g Use the small spanner provided to securely tighten all the components.
 - h Open the package containing the ion-exchange cartridges ADI. Select a cartridge and remove the plastic end caps. Place into bowl.
 - i Refit the bowl and cartridge to the de-ioniser lid. Use the large spanner to ensure a secure fit with no gap between the bowl and the lid.
 - j Now refer to OPERATING INSTRUCTIONS FOR AQUATRON DE-IONISER.

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Aquatron Filter AFH & Filter Elements AFI

Assembly Instructions

Your Aquatron Filter is a highly versatile unit which can be assembled and operated in 3 modes.

- 1 BENCH MOUNTED using the base support ring provided in the standard kit.
- 2 WALL MOUNTED using the optional wall bracket catalogue number AFD.
- 3 IN SERIES with other Aquatron de-ionisers and filters.

The assembly procedure for the bench mounted (Figure 3) and wall mounted (Figure 4) modes is almost identical to the assembly of the Aquatron De-ioniser.

Please follow the appropriate De-ioniser assembly instructions but note the following differences:

- a The bowl of the filter unit is coloured blue.
- b Plumbing components ADH/3 are used for the outlet in place of ADH/1.
- c The filter elements AFI have a paper wrapping which should be removed prior to placing in bowl.
- d Ensure the filter element is positioned centrally inside the bowl.

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
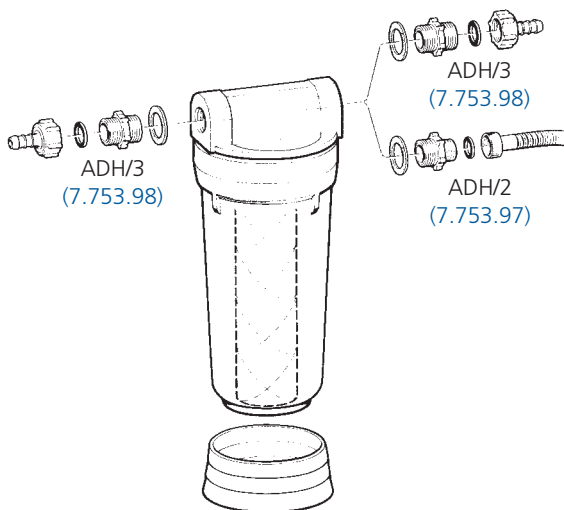
Figure 4: Filter Wall Mounted

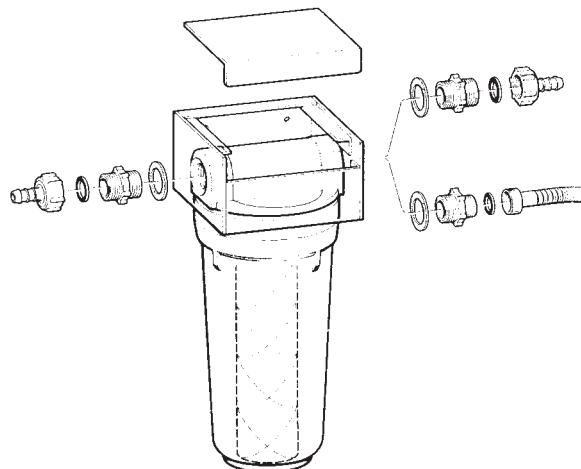Figure 4: Fixation murale du filtre Aquatron

Connecting De-ionisers and Filters in Series

It is possible to link together any number of Aquatron De-ioniser or Filter units in series. This merely requires the use of the optional plastic coupling, catalogue number ALC. One coupling is required to connect 2 units, two couplings for 3 units etc.

The following assembly procedure assumes that the series of units will be wall mounted - using the optional wall brackets AFD.

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- a Choose a suitable location for the units which provides
- a sound wall for fixing the wall brackets.
 - convenient mains water supply.
 - sufficient space to allow cartridge/filter replacement
- NB: The series of units should progress from RIGHT to LEFT. ie - The first unit accepting the mains water supply should be located on the right. The final unit in the chain will hence be on the left.

- b Take the final unit in the chain and mount on the wall by following the assembly instructions for De-ioniser and Filters on page 5. Note however the following changes:

Into the "IN" port of the de-ioniser/filter lid fit the coupling adapter ALC as shown in Figure 5.

- c Take the lid of the next unit to be assembled and position inside a wall bracket AFD. Into the "OUT" port fit the free end of the threaded adapter.
- d Offer this assembled lid to the previously wall mounted unit. Connect the $\frac{3}{4}$ BSP adapter to the $\frac{3}{4}$ BSP nut on the coupling ALC. This locates the position for the fixing of the second bracket to the wall.
- e Place the appropriate cartridges or filters into the outer bowls and fit to the CORRECT lids. Use the large spanner to ensure a secure fit with no gap between the bowl and lid.

NB: It is advisable to lubricate the rubber O ring on the lid with silicone grease. This permits easier assembly and disassembly.

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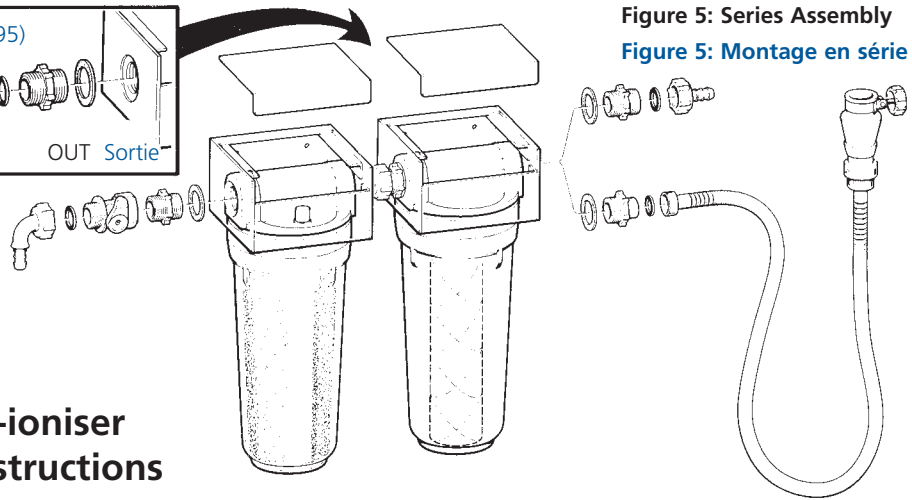
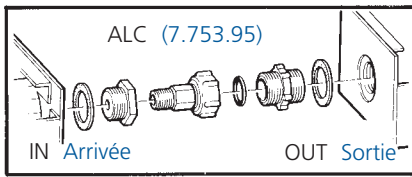
Figure 5: Series Assembly
Figure 5: Montage en série

Aquatron De-ioniser Operating Instructions

Before operating your Aquatron De-ioniser please check that:

- the maximum water temperature will not exceed 40°C.
 - the maximum water pressure will not exceed 100 psi ($7 \times 10^5 \text{ Nm}^{-2}$).
 - the plumbing fittings including hoses, will withstand the maximum pressure.
 - the pressure relief valve is CLOSED.
 - the water flow valve is OPEN.
 - a container is placed below the outlet nozzle to collect the de-ionised water.
- a Turn on the mains water supply to the de-ioniser. Observe that the unit fills with water and initially expels air from the outlet nozzle. Eventually de-ionised water emerges.
 - b Discard the first volumes of de-ionised water which is likely to be cloudy in appearance due to the presence of dissolved air.

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- c When the de-ionised water becomes clear in appearance close the water flow valve.
- d The de-ioniser is now ready for use and will produce de-ionised water on demand by opening the water flow valve.

NB: It is recommended that the de-ioniser is left under pressure (up to 100 psi) to ensure maximum cartridge efficiency.

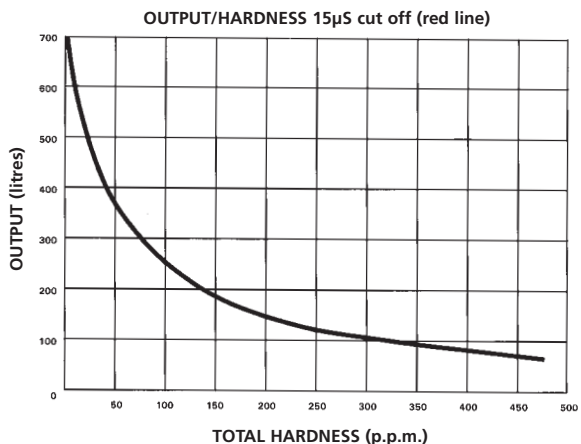
Cartridge Exhaustion/Replacement

After a period of operation the ion exchange cartridge will gradually become exhausted and the resin colour will change from green to blue.

When the colour change wave reaches the horizontal line marked on the de-ioniser bowl the cartridge should be replaced. See instructions detailed on bowl.

Notes: The quality of de-ionised water from a fresh cartridge is in the region of $1\mu\text{S}/\text{cm}$. This deteriorates to $15\mu\text{S}/\text{cm}$ when the colour change reaches the red horizontal line.

The volume of de-ionised water obtained from a cartridge is wholly dependent on the quality of the input water. Typical values are shown below:

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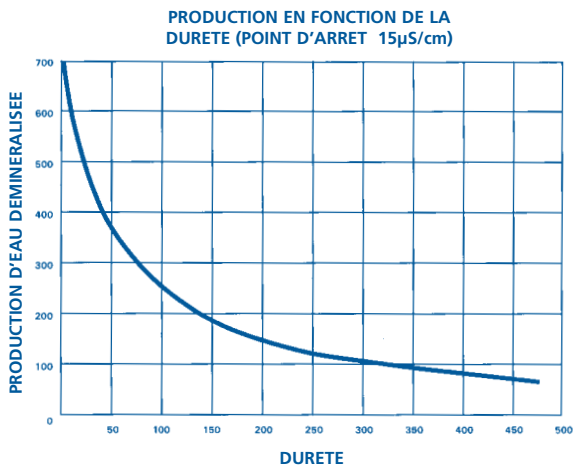
Aquatron Filter Operating Instructions

Before operating your Aquatron Filter please check that:

- the maximum water temperature will not exceed 40°C.
- the maximum water pressure will not exceed 100 psi ($7 \times 10^5 \text{ Nm}^{-2}$).
- the plumbing fittings, including hoses, will withstand the maximum pressure.
- the pressure relief valve is CLOSED.
- a container is placed below the outlet nozzle to collect the filtered water.

For operation, simply turn on the mains water supply. The filter unit will gradually fill and filtered water will emerge at the outlet nozzle.

Filter Element Replacement

After a period of operation the filter element will gradually become blocked. This is denoted by a reduction in the flow rate of filtered water.

The element should be replaced when the flowrate falls below an acceptable level and certainly BEFORE it is totally blocked.

The life of the element is wholly dependent on the quality of the input water. Experience will indicate the frequency of element replacement and once this is established it is recommended that the element be changed on a routine maintenance basis, eg. - each week, month etc.

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Accessories & Spares

Cat. No. Description

ADI	Pack of 3 ion charge cartridges.
AFI	Pack of 3 filter elements.
ADH/1	Pack of plumbing components with water flow valve.
ADH/2	Pack of plumbing components with flexible pipe and tap adapter.
ADH/3	Pack of plumbing components with plain olive connection.
AFD	Wall bracket.
ALC	Plastic coupling for mounting units in series.

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Thank you for reading this data sheet.

For pricing or for further information, please contact us at our UK Office, using the details below.

UK Office

Keison Products,

P.O. Box 2124, Chelmsford, Essex, CM1 3UP, England.

Tel: +44 (0)330 088 0560

Fax: +44 (0)1245 808399

Email: sales@keison.co.uk

Please note - Product designs and specifications are subject to change without notice. The user is responsible for determining the suitability of this product.