RB-5A, RB-12A
RB-22A
Refrigerated Baths
OPERATOR'S MANUAL

Issue 12
Date of issue:
22nd March 2011
Techne is a trademark
© Techne, 2012

CONTENTS

Refrigerated Baths User's Manual	
	page
SAFETY AND INSTALLATION	
English	4
Français	5
Deutsch	6
Español	7
INTRODUCTION	8
Operator Safety	8
Specification	9
Siting Considerations	10
Setting up the unit	10
Cleaning your unit	12
Packing	12
OPERATION	13
Precautionary Measures	13
Operating your unit	13
Daily Maintenance	13
Fault Finding	14
Service	15
ADDITIONAL INFORMATION	16
General Advice	16
General Fault finding	16
Accessories	16

Introduction																		
Please read all the information in this booklet before using the unit.																		
Warning																		
<p>HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material. Techne have taken great care in the design of these units to protect operators from hazards, but users should pay attention to the following points:</p> <ul style="list-style-type: none"> • USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS; • DO NOT put hot objects on or near combustible objects; • DO NOT operate the unit close to inflammable liquids or gases; • DO NOT place any liquid directly in your unit; • At all times USE COMMON SENSE. 																		
Operator Safety																		
All users of Techne equipment must have available the relevant literature needed to ensure their safety.																		
It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the user may be impaired.																		
All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. On some models, the cutout is adjustable and should be set to suit the application. On all other models the cutout is preset to protect the unit.																		
If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.																		
Installation																		
<ol style="list-style-type: none"> 1. All Techne units are supplied with a power cable. This may be integral or plug-in. 2. Before connecting the mains supply, check the voltage against the rating plate. Connect the mains cable to a suitable plug according to the table below. Note that the unit must be earthed to ensure proper electrical safety. <table border="0" style="margin-left: 40px;"> <tr> <td style="padding-right: 20px;"><i>Connections</i></td> <td style="padding-right: 20px;"><i>220/240V</i></td> <td><i>110/120V</i></td> </tr> <tr> <td>Live</td> <td>Brown</td> <td>Black</td> </tr> <tr> <td>Neutral</td> <td>Blue</td> <td>White</td> </tr> <tr> <td>Earth</td> <td>Green/yellow</td> <td>Green</td> </tr> </table> <p>The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable:</p> <table border="0" style="margin-left: 40px;"> <tr> <td>230V UK</td> <td>13 AMP</td> </tr> </table> <p>The fuse in the unit protects the unit and the operator.</p> <p>Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%. The rating plate is on the rear of the unit. 3. Plug the mains cable into the socket on the rear of the unit. 4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction. 5. Symbols on or near the power switch of the unit have the following meanings: <table border="0" style="margin-left: 40px;"> <tr> <td>I</td> <td>: mains switch On</td> </tr> <tr> <td>O</td> <td>: mains switch Off</td> </tr> </table> </p>	<i>Connections</i>	<i>220/240V</i>	<i>110/120V</i>	Live	Brown	Black	Neutral	Blue	White	Earth	Green/yellow	Green	230V UK	13 AMP	I	: mains switch On	O	: mains switch Off
<i>Connections</i>	<i>220/240V</i>	<i>110/120V</i>																
Live	Brown	Black																
Neutral	Blue	White																
Earth	Green/yellow	Green																
230V UK	13 AMP																	
I	: mains switch On																	
O	: mains switch Off																	
After use																		
When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.																		

Guarantee
The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.
Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.
This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.
User maintenance
NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE SIDE, FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.
THERE ARE NO USER MAINTAINABLE PARTS WITHIN THE EQUIPMENT.
In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.
<ol style="list-style-type: none"> 1. Cleaning Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool below 50° C. Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners. 2. Fuses Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.
Contact Information
For technical, sales or servicing information, contact your local Techne dealer or,
Barloworld Scientific Ltd Beacon Road, Stone Staffordshire ST15 0SA, United Kingdom Telephone: +44(0)1785 812121 Fax: +44(0)1785 813748 e-mail: equipment@barloworld-scientific.com
or,
Techne Inc, 3 Terri Lane, Suite 10, Burlington, New Jersey 08016, USA. Telephone: 609-589-2560 Toll free: 800-225-9243 ext 306 Fax: 609-589-2571 e-mail: labproducts@techneusa.com Web site: www.techneusa.com

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INTRODUCTION

Read the whole of this book before commencing work with the unit.

The book applies to a range of Techne refrigerated baths. Hence paragraphs or sections without a heading or reference apply to all units while paragraphs or sections with headings apply to specific units. For other refrigerated units there is a separate booklet.

The Techne fridge units are designed to cool liquids to temperatures around ambient and below by extracting heat continuously. The temperature is controlled by a temperature controller with a heater adding heat as necessary. Depending upon the circuit and the application this could be a Techne Thermoregulator, TE-10 series or TU-20 series. Each unit, fridge or controller, has its particular application which should be considered when setting up a temperature controlled system.

The refrigerated baths have integral baths and have refrigeration coils within the bath. They are designed to work with Techne thermoregulators. The Fridge Bath works best when there is a good flow of liquid over the coil.

Operator Safety

All users of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures.

Techne refrigerated units have been designed to conform to international safety requirements.

If there is any doubt relating to the proper use of this equipment, the staff at Techne will be pleased to assist you.

Specification

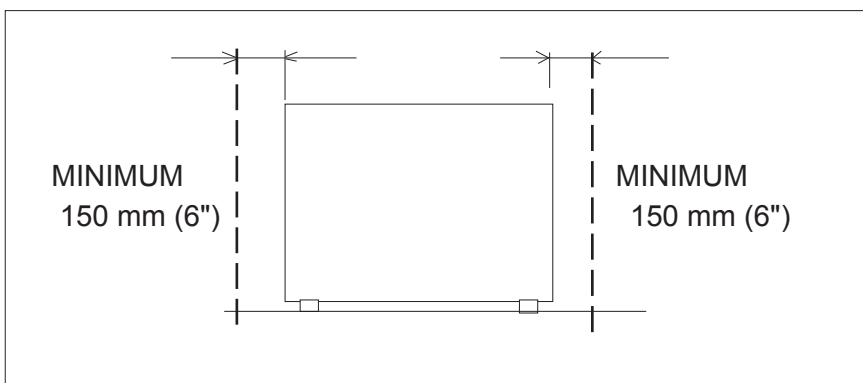
The minimum temperatures quoted in the specification are achieved in an ambient of 20 °C with the appropriate liquid in the integral bath. However many things will affect the minimum obtainable temperature, such as: ambient temperature, insulation and length of circulation system, condition of the working liquid, heat input and so on.

	RB-5A	RB-12A	RB-22A
Minimum temperature	-20 °C	-35°C	-30°C
Cooling capacity at:			
20 °C	145 W	240W	240W
0 °C	145 W	240W	240W
-10 °C	110 W	230W	230W
Nominal dimensions and weights			
Overall:			
Length	430 mm	430 mm	430 mm
Width	250 mm	370 mm	370 mm
Height	566 mm	610 mm	690 mm
Bath opening:			
Length	192 mm	208 mm	208 mm
Width	151 mm	300 mm	300 mm
Depth	200 mm	150 mm	225 mm
Maximum working capacity	7 litres	11.6 litres	22 litres
Net Weight	23 kg	44 kg	49 kg

* Including the Thermoregulator

Siting Considerations

All the units are air cooled, air being pulled in through the front and expelled from the rear. The ventilation grilles should never be obstructed, therefore, as this will detract from the performance. Also the fridge units should never be situated in an enclosed space or near an extractor fan, this too will detract from the performance.

Setting up the unit

1. Check the voltage marked on the unit's rating plate against that of the mains supply. Connect the mains cable to a suitable plug as in the following table. **Note that the unit must be earthed to ensure proper electrical safety.**

<i>Connections</i>	<i>230V</i>	<i>110/120V</i>
Live	brown	black
Neutral	blue	white
Earth	yellow and green	green

The fused plug supplied on the mains cable for use in the UK with the RB-5A is fitted with the following fuse to protect the cable.

<i>Supply</i>	<i>fuse fitted</i>
240V	5 Amp

For the RB-12A and RB-22A, if a fused plug is fitted then a fuse of the following value must be fitted

<i>Supply</i>	<i>fuse fitted</i>
240V	13 Amp

The fuse in the unit protects the unit and the operator.

Note that units marked 230V on the rating plate work equally well at 220V; units marked 120V on the rating plate work equally well at 110V.

2 Recommended liquids:

When working at low temperatures always use a liquid suitable for the lowest temperature at which the unit is capable of achieving, see the list of recommended liquids below. When using a flow cooler up to 30°C in an enclosed circuit use the recommended liquid.

For the **RB-12A** and **RB-22A** units:


<i>Temperature</i>	<i>Liquid</i>
-35 °C to 0 °C *	40% water 40% anti-freeze 20% alcohol

For the **RB-5A** units:

<i>Temperature</i>	<i>Liquid</i>
-20 °C to 30 °C *	50% water 50% anti-freeze.

* **Warning:** Above 0°C alcohol may evaporate from the liquid giving a health and safety hazard and also deteriorating the mix. Check with health and safety procedures before using above 0°C. Above 30°C anti-freeze may give off dangerous fumes. Again check with health and safety procedures before using above 30°C. There is no liquid which will cover the full range of working temperatures. The RB-12A & RB-22A units are designed to work up to 40°C with occasional peaks to 100°C. The RB-5A can be used up to 150°C but unit should be turned off when operating at temperatures above 80°C.

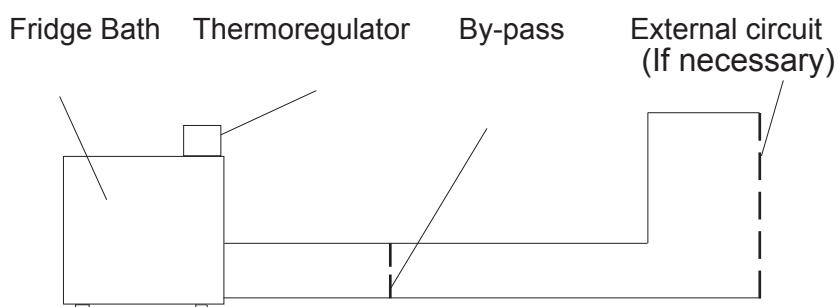- Hoses should be as short as possible and secured with clips.
- The whole system should be insulated as much as possible.
- Set up the bath on a flat surface.
- For fixing the thermoregulator to the bridge, see the thermoregulator manual.
- If external circulation has been chosen suitable hoses should be fitted to the outlet and return pipes of the thermoregulator. Always securely clip the hoses in place. A suitable hose is one of 16mm bore and capable of withstanding both the temperature of operation and the liquid being used. 12mm bore hose may be used if the hose is soft such as silicone rubber.
- If the system is of a small bore or the viscosity of the liquid prevents the flow being maintained, a by-pass should be incorporated in the system.

- 9 Fill the bath to between the minimum and maximum levels which are 65mm and 20mm from the top of the bath.
- 10 Fridge baths are supplied with a lid and this should be used whenever possible..

Cleaning your unit

Before cleaning your unit, disconnect from the power supply and allow to cool to ambient temperature.

You can clean the case of the refrigerated unit with a cloth dipped in water or ethanol (methanol or formaldehyde can also be used). No part of the case or cover should be immersed in the solvents.

Do not use acetone or abrasive cleaners.

Packing

When you receive your new refrigerated bath, remove all the packing. Check the unit for any transit damage. If there is any transit damage at all inform Techne Sales Department or your Techne dealer immediately.

When unpacking the unit, check that the following have been removed from the packing: a power cable (RB-5A only); a decontamination certificate and a Guarantee Card.

Check that the unit is the correct rating for your power supply by looking on the Rating Plate which is on the rear of the unit.

Within the guarantee period, shown on the Guarantee Card, we undertake to supply replacements free of charge for parts which may, on examination, prove to be defective, provided that the defect is not the result of misuse, accident or negligence. Any instrument requiring service under this guarantee should be sent to the supplier through whom it was purchased, or, in the case of difficulty, it should be carefully packed in its original packing and consigned, carriage paid, to us. Techne takes no responsibility for returned goods damaged in transit.

Returned goods will not be processed without a Returns Authorisation Number. Call + (44) (0)1785 810475 for a number. On all correspondence, please quote the Serial Number in full and/or the Sales Order Number. **Please write the Returns Number on the outside of any packing.**

Precautionary Measures

- 1 Do not operate any of the fridge units without liquid in the circuit. Ensure that the evaporation coil is covered.
- 2 Freezing may occur if there is insufficient circulation or the incorrect liquid is used. While this will not harm the unit some time may pass before the unit can be used again. For this reason there should always be circulation in or around the coil either in the bath. Loss of performance will occur if freezing takes place.
- 3 To avoid thermal shock do not fill a cold bath with, liquids above 40°C.
- 4 Although the practice is not recommended fridge units may occasionally be subjected to temperatures up to 100°C. This is provided that they are only being used thus intermittently, for example for cooling down to temperatures below 40°C.
- 5 IN NO CIRCUMSTANCES MAY A REFRIGERATED UNIT BE SUBJECTED TO BATH/LIQUID TEMPERATURES ABOVE 100°C.

Operating your unit

- 1 Switch on the thermoregulator as instructed in the appropriate book so that good circulation is established.
- 2 Switch the fridge unit on by pressing the power switch. The switch and the indicator on the front will light up. The fan and the compressor will start. It will take a little while for cooling to begin.
- 3 Remember that as the Fridge units will only extract heat, a controller, such as a TE-10 or TU-20 thermoregulator, will need to be incorporated in the system. Temperatures are set on and controlled by the temperature controller, see the appropriate book.

Daily Maintenance

As mentioned earlier, it is imperative that the air flow is not impeded in any way. It is important that the grilles are kept clean and clear at all times. Even dust and fluff can impair the performance, therefore the units must be used in a clean environment or they should be regularly cleaned.

Fault Finding

- 1 Check that the airways have not been restricted either by something placed in front of them or through build up of dust and fluff on the grilles, etc.
- 2 Loss of performance is most often caused by deterioration of the working liquid. As a first check always mix up a new batch of liquid and re-test the unit. NEVER top up the mix with water always use a quantity of the proper mix.
- 3 The next most common cause of poor performance is that the liquid is not being circulated over the coil or through the heat exchanger properly. First check that the pump unit is not blocked either by foreign matter or by balls from a ball blanket or in any other way. Also check that the pipe system, if you are using the fridge unit for an external system, is not kinked or restricted in any way.
- 4 If the liquid or the circulation is not the cause of the poor performance then there may be a number of other causes.
- 5 Check that the internal fan is running by feeling if there is air blowing when the unit is on. If not then the unit needs servicing as below.
- 6 Check that the compressor is running by listening for it to start and stop as the unit gets to its lowest temperature. If not then the unit needs servicing as below.
- 7 Last of all the unit may have lost gas and needs servicing as below
- 8 For detailed repair instructions of the thermoregulator refer to the appropriate model instructions.

NOTE: Before any work is undertaken on the thermoregulator it MUST be removed from the bath.

Service

- 1 Complete servicing and repair of Techne products is available at Techne see the contact information on pages 3 or 4.
- 2 A Service Manual is available and servicing may be undertaken by any competent refrigeration engineer.

The following table gives a guide as to the gas charge for each of the units:

<i>Unit</i>	<i>Refrigerant §</i>	<i>Charge *</i>
RB-5A	R134a	85 gm
RB-12A	R404a	250 gm
RB-22A	R404a	250 gm

- * There have been several improvements in design which have varied the charge. To increase the cool-down rate increase the charge slightly; to reach a lower temperature reduce the charge slightly. Adjust the charge during testing.
- § Early units were charged with R12 or R502 please consult Techne for advice on repair of older units. Units charged with R404a or R134a have GN after the Serial Number, eg 12345-13gn.

ADDITIONAL INFORMATION

Brief fault finding notes and lists of accessories are given in this section.

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL.

REMOVING THE OUTER COVERS OR BASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES AND A FAN.

THERE ARE NO USER SERVICEABLE PARTS WITHIN THE EQUIPMENT.

General advice

In the unlikely event that you experience any problems with your refrigerated unit which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier.

Fuses

If the power light on the power switch is lit, the fuse may have blown. Make sure there is no external cause (such as a faulty plug or lead).

Fuses should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.


Replacement fuse numbers are as follows, and can be ordered from Techne or thier agents;


Number	Rating		Unit voltage
RB-5A			
6008444	T4A	4 amp time delay	230V
6501239	T8A	8 amp time delay	120V
RB-12A or RB-22A			
6008442	T8A	4 amp time delay	230V
6008443	T12.5A	8 amp time delay	120V

Accessories

There are many accessories, these are constantly being updated and details can be found in Techne Sales literature. They include:

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- Temperature Program Software;
- Cooling Water Pack;
- Refrigeration Control Pack;
- Test Tube Racks;
- and so on.

Thank you for reading this data sheet.

For pricing or for further information, please contact us at our UK Office, using the details below.

UK Office

Keison Products,

P.O. Box 2124, Chelmsford, Essex, CM1 3UP, England.

Tel: +44 (0)330 088 0560

Fax: +44 (0)1245 808399

Email: sales@keison.co.uk

Please note - Product designs and specifications are subject to change without notice. The user is responsible for determining the suitability of this product.