
CH-5 and PC-5
Chiller unit and Pump
OPERATOR'S MANUAL

Issue 12.
Date of issue:
3rd May 2005
Techne is a trademark
© Techne , 2003

CONTENTS**CH-5 Chiller User's Manual**

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	page
SAFETY AND INSTALLATION	
English	4
Français	5
Deutsch	6
Español	7
INTRODUCTION	8
Specification	9
Siting Considerations	10
Setting up the unit	10
OPERATION	12
Precautionary Measures	12
Operating your unit	13
Setting the temperature	13
Running the unit	13
ADDITIONAL INFORMATION	14
Spare Parts	16
Accessories	16

NOTES

Introduction

Please read all the information in this booklet before using the unit.

Warning

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material. Techne have taken great care in the design of these units to protect operators from hazards, but users should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS;
- DO NOT put hot objects on or near combustible objects;
- DO NOT operate the unit close to inflammable liquids or gases;
- DO NOT place any liquid directly in your unit;
- At all times USE COMMON SENSE.

Operator Safety

All users of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the user may be impaired.

All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. On some models, the cutout is adjustable and should be set to suit the application. On all other models the cutout is preset to protect the unit.

If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.

Installation

1. All Techne units are supplied with a power cable. This may be integral or plug-in.
2. Before connecting the mains supply, check the voltage against the rating plate. Connect the mains cable to a suitable plug according to the table below. **Note that the unit must be earthed to ensure proper electrical safety.**

Connections	220/240V	110/120V
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable:

230V UK	13 AMP
---------	--------

The fuse in the unit protects the unit and the operator. Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%. The rating plate is on the rear of the unit.

3. Plug the mains cable into the socket on the rear of the unit.
4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
5. Symbols on or near the power switch of the unit have the following meanings:

I	: mains switch On
O	: mains switch Off

After use

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.

Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.

Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.

This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.

User maintenance

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL.

REMOVING THE SIDE, FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.

THERE ARE NO USER MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

1. **Cleaning**

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool below 50° C. Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

2. **Fuses**

Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

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
INTRODUCTION

Read the whole of this book before commencing work with the unit.

The Techne CH-5 fridge units are designed to work with the Techne PC-5 at temperatures between +4°C and +15°C using demineralised water. Each unit is fitted with a thermostat which controls the temperature of liquid in the bath. The bath has a refrigeration coil which extracts heat from the liquid. They work most efficiently when there is a good flow of liquid over the coil.

The PC-5 is specifically designed to work with the CH-5. It has a double outlet to deliver chilled water and a connection to return the water to the bath. Care must be taken when installing the units.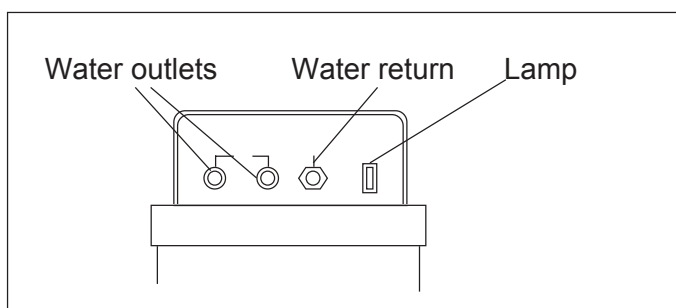

Specification

The minimum temperatures quoted in the specification are achieved in an ambient of 20°C with the appropriate liquid in the integral bath. However many things will affect the minimum obtainable temperature, such as: ambient temperature, insulation and length of circulation system, condition of the working liquid, heat input and so on.

CH-5

Minimum temperature achievable	+4°C
Cooling capacity at: 20°C	400 W
4°C	400 W

Nominal dimensions and weights

Overall:

Length	430 mm
Width	235 mm
Height	402 mm

Bath Opening:

Length	300 mm
Width	150 mm
Depth	150 mm

Net Weight	23 kg
------------	-------

PC-5

Pump Capacity	20 l/min
Pump Head	6 meters of water


Nominal dimensions and weights

Overall:

Length	380 mm
Width	235 mm
Height above bath	122 mm
Extension into bath	80 mm
Net Weight	3.5 Kg

Siting Considerations

The units are air cooled, air being pulled in through the front and expelled from the rear. The ventilation grilles should never be obstructed, therefore, as this will detract from the performance. Also the fridge units should never be situated in an enclosed space or near an extractor fan, this too will detract from the performance.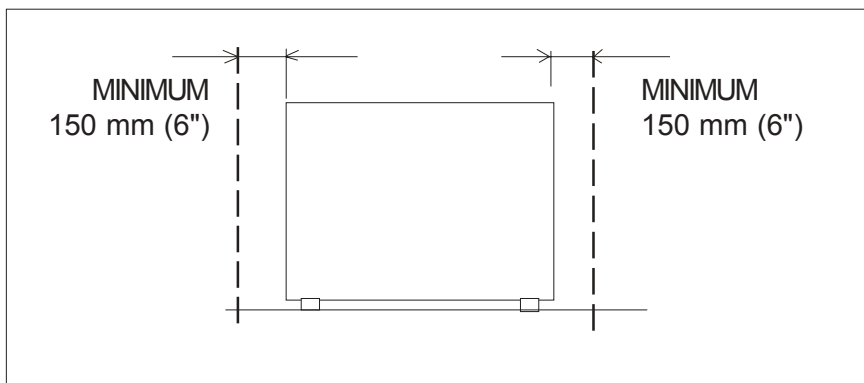

Setting up the unit

1. Check the voltage marked on the unit's rating plate against that of the mains supply. Connect the mains cable to a suitable plug as in the following table. **Note that the unit must be earthed to ensure proper electrical safety.**

<i>Connections</i>	<i>220/240V</i>	<i>110/120V</i>
Live	brown	black
Neutral	blue	white
Earth	yellow and green green	


The fused plug supplied on the mains cable for use in the UK with the RB-5A is fitted with the following fuse to protect the cable.

<i>Supply</i>	<i>fuse fitted</i>
240V	5 Amp


The fuse in the unit protects the unit and the operator.

Note that units marked 240V on the rating plate work equally well at 220V; units marked 120V on the rating plate work equally well at 110V.

- 2 Set up the chiller on a level surface. Use demineralised water to fill the bath and cover the cooling coil. Demineralised water is preferred to reduce the formation of scale. If scale should form, use only mild de-scaling agents to remove it. DO NOT attempt to hammer, chip or scrape the deposits from the surface of the bath or the coil.
- 3 Place the PC-5 on the CH-5 and position with the studs on the chassis of the pump unit.
- 4 Connect the PC-5 to the unit you wish to chill. Ensure that all connections are secure and will withstand the pressure from the pump.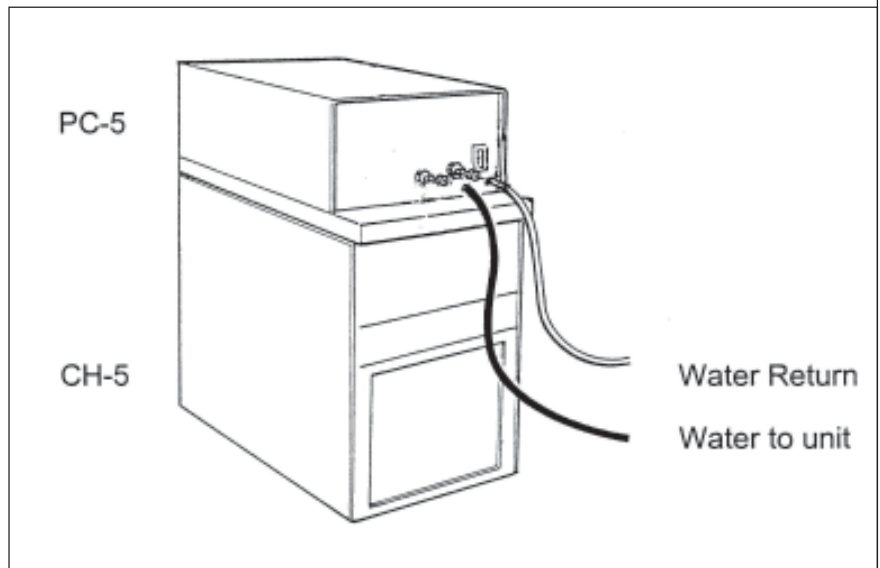

- 5 Plug the in line plug on the PC-5 into the socket on the back of the CH-5.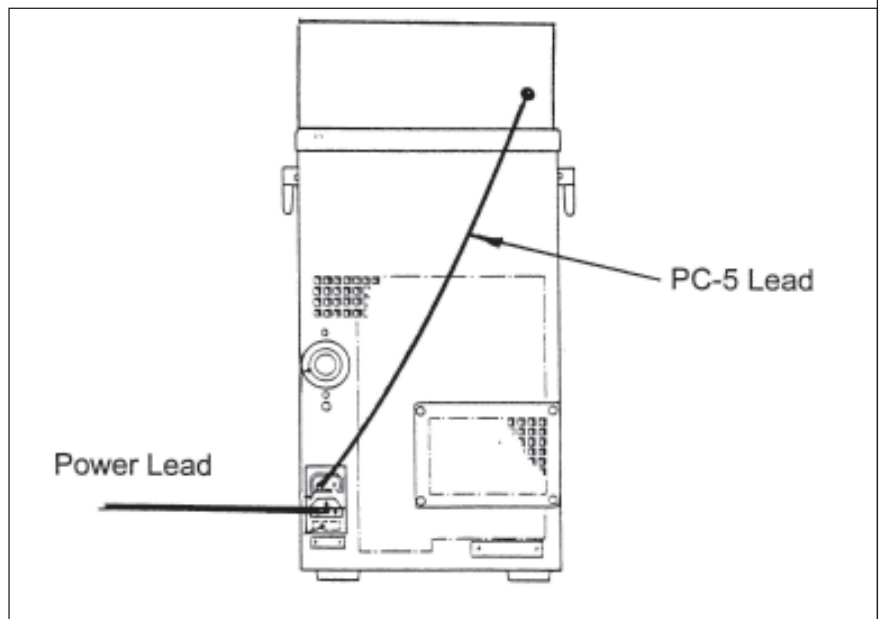

- 6 Insulate all the connection pipes to minimise condensation when working at low temperatures.
- 7 Both the PC-5 and the CH-5 are now ready to run.

OPERATION


Precautionary Measures

- 1 Do not operate the fridge unit without liquid in the bath such that the cooling coil is covered.
- 2 Do not operate the fridge unit without the pump unit working to give good circulation in the bath as freezing up may occur. While this will not harm the unit some time may pass before the unit can be used again. Loss of performance will occur if freezing does take place.
- 3 Do not operate the pump without proper connections both out of the bath and returning to the bath.
- 4 Although the practice is not recommended fridge units may occasionally be subjected to temperatures up to 100°C. This is provided that they are only being used thus intermittently, for example for cooling down to temperatures below 40°C.
- 5 IN NO CIRCUMSTANCES MAY A REFRIGERATED UNIT BE SUBJECTED TO BATH/LIQUID TEMPERATURES ABOVE 100°C.
- 6 IF YOUR UNIT SHOULD ICE UP DO NOT TRY TO FORCE THE PUMP UNIT FROM THE BATH THIS MAY DAMAGE BOTH THE CHILLER AND THE PUMP. Allow the bath to de-freeze naturally and then lift the pump from the water. Inspect for damage then adjust the unit as under Running the Unit.
- 7 NEVER run the chiller unit without the pump running as this may result in localised freezing.

Operating Your Unit

Setting the temperature.

The temperature of the recirculating liquid is set by the thermostat dial on the rear of the unit.

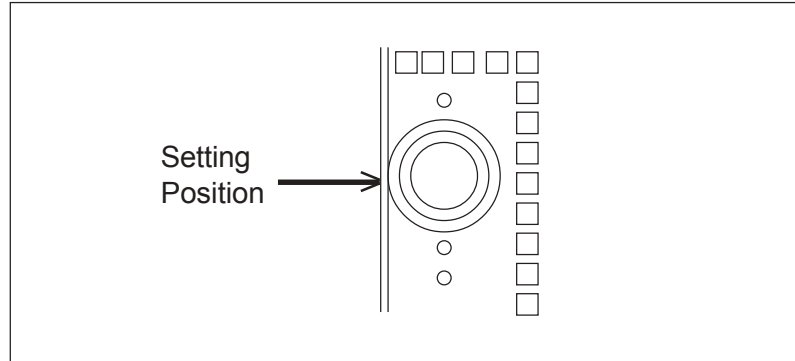Setting the dial to 1, gives a temperature of approximately 15°C.

Setting the dial to 3, gives a temperature of approximately +4°C.

If there is no flow it is quite likely that the water has frozen. Do NOT attempt to lift the pump from the bath as this can damage the coil. Wait for the ice to turn to water. There is no need to remove the pump except to clean the bath and refill it with water. The water will freeze if, you set the dial to a higher number than 4, and/or, if there is no heat for the water to extract from the unit being feed with chilled water. The CH-5 is a robust unit if you take these precautions.

Running the Unit

- 1 Set the thermostat dial to 3 or 4.
- 2 Switch the Chiller unit on by pressing the power switch. The switch and the indicators on the front of the CH-5 and PC-5 will illuminate. The fan and the compressor will start in the Chiller and the pump will start in the Pump unit. It will take a little while for cooling to begin.
- 3 The Pump will run continuously.
- 4 Measure the temperature of the water and adjust the dial until the desired temperature is attained. Remember that if the water freezes, do not force the pump from the bath.

ADDITIONAL INFORMATION

Brief fault finding notes and lists of accessories are given in this section.

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. **REMOVING THE OUTER COVERS OR BASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.**

THERE ARE NO USER SERVICEABLE PARTS WITHIN THE EQUIPMENT.

General advice

In the unlikely event that you experience any problems with your refrigerated unit which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier.

Packing

When you receive your unit, make sure you keep the original packing in case you ever need to return it for service or repair.

The unit must be transported in the original packing to avoid damage. Techne accepts no responsibility for damage incurred unless the unit is correctly packed and transported in this way. General fault finding

Fuses

If the power light on the power switch is lit, the fuse may have blown. Make sure there is no external cause (such as a faulty plug or lead).

Fuses should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

Cleaning your unit

Before cleaning your unit, disconnect from the power supply and allow to cool to ambient temperature.

You can clean the case of the refrigerated unit with a cloth dipped in water or ethanol (methanol or formaldehyde can also be used). No part of the case or cover should be immersed in the solvents.

Do not use acetone or abrasive cleaners.

Daily Maintenance

As mentioned earlier, it is imperative that the air flow is not impeded in any way. It is important that the grilles are kept clean and clear at all times. Even dust and fluff can impair the performance, therefore the units must be used in a clean environment or they should be regularly cleaned.

Fault Finding

- 1 Check that the airways have not been restricted either by something placed in front of them or through build up of dust and fluff on the grilles, etc.
- 2 The most common cause of poor performance is that the liquid is not being circulated over the coil. First check that the pump unit is not blocked either by foreign matter or by balls from a ball blanket or in any other way. Also check that the pipe system, if you are using the fridge unit for an external system, is not kinked or restricted in any way.
- 4 If the liquid or the circulation is not the cause of the poor performance then there may be a number of other causes.
- 5 Check that the internal fan is running by feeling if there is air blowing when the unit is on. If not then the unit needs servicing as below.
- 6 Check that the compressor is running by listening for it to start and stop as the unit gets to its lowest temperature. If not then the unit needs servicing as below.
- 7 Last of all the unit may have lost gas and needs servicing as below

Service

- 1 Complete servicing and repair of Techne products is available at Techne see the contact information on page 4.
- 2 A Service Manual is available and servicing may be undertaken by any competent refrigeration engineer.

The following table gives a guide as to the gas charge for each of the units:

<i>Unit</i>	<i>Refrigerant §</i>	<i>Charge *</i>
CH-5	R134a	95 gm

- * There have been several improvements in design which have varied the charge. To increase the cool-down rate increase the charge slightly; to reach a lower temperature reduce the charge slightly. Adjust the charge during testing.
- § Early units were charged with R12 please consult Techne for advice on repair of older units. Units charged with R134a have GN after the Serial Number, eg 12345/5 GN.

Accessories

There are many accessories, these are constantly being updated and details can be found in Techne Sales literature. They include:

- Boost Units;
- Temperature Programmers;
- Cooling Water Pack;
- Test Tube Racks and so on.

Thank you for reading this data sheet.

For pricing or for further information, please contact us at our UK Office, using the details below.

UK Office

Keison Products,

P.O. Box 2124, Chelmsford, Essex, CM1 3UP, England.

Tel: +44 (0)330 088 0560

Fax: +44 (0)1245 808399

Email: sales@keison.co.uk

Please note - Product designs and specifications are subject to change without notice. The user is responsible for determining the suitability of this product.